

DH6 / DH6-LC Cleaning of Cooling Components

Rev. 1.1 - 06. Mar. 2026

This Service Note explains the process of cleaning the cooling components of a DH6 or DH6-LC if needed.

Safety Instructions

- Remove any Ozone or overpressure from the tubing
- Disconnect the device from mains power
- Attach your anti-static wrist strap to ground, only – if this is not possible: at least eliminate any electrostatic charge by touching a grounded surface before opening the device
- Make sure to wear eye protection




Description

1. Open up the instrument's back panel with a 2mm Allen Wrench. (6 screws, on older version 4 screws)



2. Unplug the fan connector from the fan.

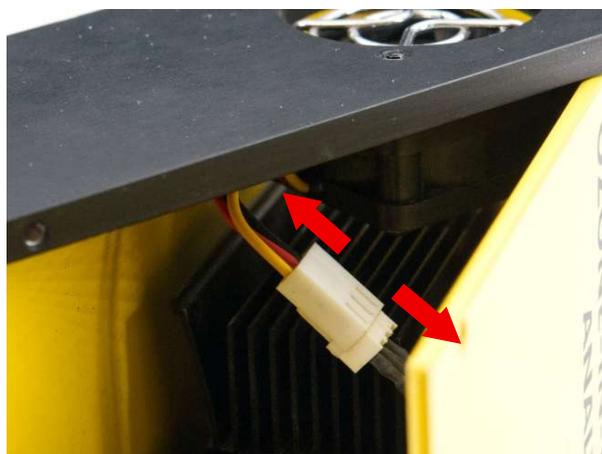

3. The fan and the cooling fins can be cleaned with lint-free cloth and compressed air. Use only low-pressure airflow and avoid excessive air volume, as this may damage the fan. Make sure to wear eye protection.



4. Repeat the steps in reverse order, to put the DH6 / DH6-LC back together.
Do not forget to reattach the fan connector.

If any problems occur, do not hesitate to contact us:

BMT Messtechnik GmbH, Hamburger Str. 19, D-14532 Stahnsdorf, Germany
Tel.: +49-3329-696 77-0
service@bmt-berlin.de

Your BMT Service Team